2025/11/05 13:15 1/3 About

## **About**



The ThermoFisher Laboratory Automation System (LAS) serves as the cornerstone technology for the Living Biofoundry, enabling execution of automated customized synthetic biology and workflows at >500 samples-per-week. The LAS is equipped with over of over 10 functional instrumental components, including: a state-of-the-art SpinnakerTM microplate robot, automated incubators, reagent dispensers, thermal cyclers, plate sealer, and carousels/racks that are seamlessly integrated through the MOMENTUM application programming interface that is fully-compatible with laboratory information management systems (LIMS).

# **Equipment Available**

## **Thermo Scientific Multidrop Combi**



#### **About**

The Multidrop Combi is an automatic, programmable, eight-channel microplate bulk reagent dispenser for microvolume dispensing. It can dispense 1 to 8 different reagents from external liquid reservoirs into different rows and can be used in several applications, for example, in drug discovery/highthroughput screening, genomic, proteomic and cell-based assays, and ELISA. It has a volume range of 0.5 to 2500  $\mu$ l for 96-well plates of heights 5-50 mm.

#### **User Notes**

**Before Use**: Place the dispensing cassette tubes in proper position. Ensure that the waste containers are not full.

**After Use**: Flush out the tubing with deionized distilled water. Empty out waste containers, disposing of the contents as appropriate. Place the dispensing cassette tubes in rest position. For more information please read:

#### multidrop combi.pdf

## **BioTek ELx405 Microplate Washer**



#### **About**

The BioTek ELx405 is the first deep well plate washer that aspirates and dispenses into 96- and 384-well deep well and standard height microplates. It is a versatile, high-performance laboratory instrument designed for automated washing of microplates, supporting a wide range of applications such as ELISA, cell-based assays, immunoassays, etc.

Key Features:

**Plate Compatibility**: Washes 384-well and 96- well microplates in addition to 24-well and 96-well deep well plates (up to 50 mm in height)

### **Washing Technology:**

Utilizes a patented Dual-Action™ manifold, enabling independent control of dispense and aspiration for precise washing, overfill washing, and overflow protection.

Low-flow rates and 20° angled dispensing are available, making it especially suitable for gentle washing of adherent cell-based assays.

### Fluid Handling:

Internal positive displacement pump for accurate fluid delivery, supporting volumes from 50 to 3000  $\mu$ L/well.

Optional buffer switching allows for up to four different wash buffers to be used automatically in a single protocol.

#### **User Notes**

**Before Use**: Ensure that the waste containers are not full.

**After Use**: Flush out washer with deionized water. Empty out waste containers, disposing of the contents as appropriate. For more information please read:

elx405cleaning.pdf

## **Automated Thermal Cycler**

## **About**

The Automated Thermal cycler (ATC) is capable of performing PCRs (polymerase chain reactions) at a high-throughput scale. Other features include:

- 1. Automated lid critical for hands-free operation
- 2. Automated plate ejectors to ensure consistent plate positioning
- 3. Heated cover to protect the plate nest, minimizing evaporation and enabling thermal contact

2025/11/05 13:15 3/3 About

# **Reference documentation**

momentum\_6.1\_training\_module\_4\_v3\_spinnaker\_advanced.pdf

spinnaker\_user\_guide.pdf

30124664\_ifu\_spark\_ref\_english\_v2\_0.pdf

operating\_instructions\_cytomat\_44.pdf

elx405 user manual sd-xb000358.pdf

440140 - azenta xpeel user manual.pdf

multidropcombimanual.pdf

From:

https://bpm-wiki.cnsi.ucsb.edu/ - NSF BioPACIFIC MIP Wiki

Permanent link:

https://bpm-wiki.cnsi.ucsb.edu/doku.php?id=lbf-las&rev=1751069605

Last update: 2025/06/28 00:13

