

# Harrick Plasma Cleaner PDC-32G

Plasme Cleaner



**Tool Type:** "Cleaning Tool"

**Location:** "Microfluidics Lab"

| Supervisor            | Tool Lead      |
|-----------------------|----------------|
| David Bothman         | "WW Name"      |
| (805) 893-4125        | (###) ###-#### |
| bothman@cnsi.ucsb.edu | "WW Email"     |

**Description:** "TEXT HERE"

**Manufacturer:** "TEXT HERE"

## About

The Harrick PDC-32G is a compact table top plasma cleaner for nanoscale surface cleaning and activation of small samples. The attached PlasmaFlo gas mixer allows for quantitative control of up to two different process gasses and monitoring of the vacuumed pressure within the chamber.

## Training Documentation

[Plasma Cleaner SOP](#)

## Detailed Specifications

- 3 stage adjustable RF power settings
- Max RF power: 18W

- Chamber size: 3" diameter x 6.5" Pyrex chamber
  - Valve threads: 1/8 NPT
- 

## Recipes

Brian Dincau found that the following recipe does a good job enhancing PDMS-PDMS bonding:

- Power - High
- Pressure - 450 mTorr
- Duration - 30s
- Nitrogen flow -

## Safety Concerns

---

## Reference Documentation

[harrick-plasma-product-information-web.pdf](#)

[plasma\\_cleaner\\_manual\\_32g.pdf](#)

[ufl\\_harrick\\_plasma\\_operating\\_instructions.docx](#)

---

From:

<https://bpm-wiki.cnsi.ucsb.edu/> - NSF BioPACIFIC MIP Wiki

Permanent link:

[https://bpm-wiki.cnsi.ucsb.edu/doku.php?id=plasma\\_cleaner&rev=1603302507](https://bpm-wiki.cnsi.ucsb.edu/doku.php?id=plasma_cleaner&rev=1603302507)

Last update: **2020/10/21 17:48**

