

# CNSI Microfluidics Laser Cutting Data

## Material Vendors

Acrylic sheet, optically clear, from McMaster Carr: part number 8560K239  
<http://www.mcmaster.com/#catalog/119/3558/=molz4w>

## Cutting Parameters

### Acrylic Sheet

Cutting 1/8", 100% power, 0.4 ipm, 2000 Hz, (Dave Bothman, 5/5/13 - clean cut)

Cutting 1/4", 100% power, 0.15 ipm, 1000 Hz, z offset -.125, (Dave Bothman, 5/5/13 - clean cut)

### PDMS

250um PDMS with mylar backing on back side and removed from front

Cutting 42% power, 2.0 ips, 5000Hz (Peter Mage, 5/9/13 - some flash and soot - recipe needs to be refined)

### Double Sticky Tape

25% power, 3 ips

## Guideline Cutting Recipes

Heading 1	Heading 2	Heading 3
Row 1 Col 1	Row 1 Col 2	Row 1 Col 3
Row 2 Col 1	some colspan (note the double pipe)	
Row 3 Col 1	Row 3 Col 2	Row 3 Col 3

- Recipes from Vincent

Heading 1	Heading 2	Heading 3
Row 1 Col 1	Row 1 Col 2	Row 1 Col 3
Row 2 Col 1	some colspan (note the double pipe)	
Row 3 Col 1	Row 3 Col 2	Row 3 Col 3

- Recipes from Rachel

Heading 1	Heading 2	Heading 3
Row 1 Col 1	Row 1 Col 2	Row 1 Col 3
Row 2 Col 1	some colspan (note the double pipe)	
Row 3 Col 1	Row 3 Col 2	Row 3 Col 3

- Recipes from Ola

Heading 1	Heading 2	Heading 3
Row 1 Col 1	Row 1 Col 2	Row 1 Col 3
Row 2 Col 1	some colspan (note the double pipe)	
Row 3 Col 1	Row 3 Col 2	Row 3 Col 3

- Recipes from Beach
- 

From: <https://bpm-wiki.cnsi.ucsb.edu/dokuwiki/> - NSF BioPACIFIC MIP Wiki

Permanent link: [https://bpm-wiki.cnsi.ucsb.edu/dokuwiki/doku.php?id=laser\\_cutting\\_data&rev=1405545877](https://bpm-wiki.cnsi.ucsb.edu/dokuwiki/doku.php?id=laser_cutting_data&rev=1405545877)

Last update: **2014/07/16 21:24**

