
2025/10/16 00:02 1/3 SMaLL Printer

SMaLL Printer

SMaLL Printer

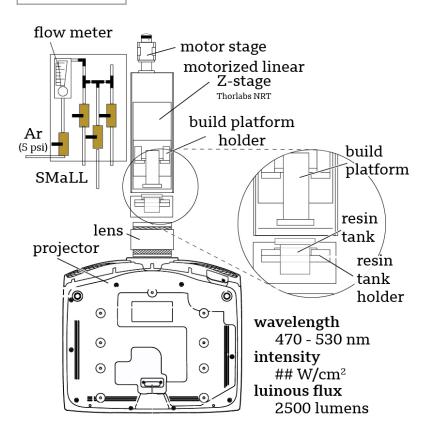

Tool Type: "3D Printer"
Location: "Elings Hall 2436"

Supervisor	Tool Lead
Juan Manuel Urueña	
jmuruena@ucsb.edu	"WW Email"

Description: "Solution Mask Liquid Lithography (SMaLL)"

Manufacturer: "In-House"

About

SMaLL is capable of creating unique structures with mechanically and chemically distinct regions. By using tunable, visible wavelength photoswitches, in combination with orthogonal crosslinking reactions, a single resin formulation can be employed for the facile, one-step printing of complex, bioinspired structures. The fabrication of soft joints and brick-and-motar architectures showcase the utility of SMaLL for printing all-polymer composites with well-defined regions for different mechanical and chemical properties.

Last update: 2021/12/13 21:26

Detailed Specifications

wavelength: 420 - 700 nmluminous flux: 2,500 lumens

Projector: Epson PowerLite Home Cinema 5040UB
1080p native resolution with 4K Enhancement

Safety Concerns

Read the manufactures manual before first use. If the BIO X Cellink acts in a way that is not described by the manual, turn off the printer and contact CELLINK.


- Never place your finger near the machine until all parts have stopped moving. Moving parts can cause serious injury
- Never clean or service the printer while it is on
- The printer uses UV light for sterilization and curing. Never look directly at UV light nor expose skin. Serious injury may result from exposure
- The printer has heated surfaces that can reach temperature up to 250 degrees Celsius. Never touch these surfaces when using the heating function. Allow things to cool before opening, touching the printer
- Always ensure that equipment is correctly mounted before use. Imporperly mounted print beds, printheads, cartridges, calbes liquid spouts and air spouts can be dangerous. If any equipment appears damaged, turn off the printer, unplug all connections and contact CELLINK.

Operating Procedures

Insert Text Here!

Reference Documentation

2025/10/16 00:02 3/3 SMaLL Printer

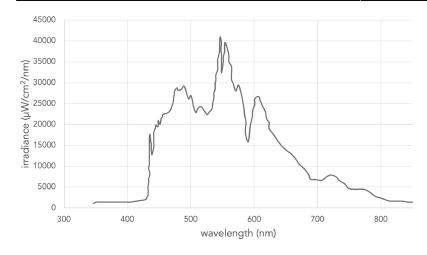

creationworkshop.pdf

small_projector_manual.pdf

Training Documentation

small_projector_manual.pdf

From

https://bpm-wiki.cnsi.ucsb.edu/dokuwiki/ - NSF BioPACIFIC MIP Wiki

Permanent link:

https://bpm-wiki.cnsi.ucsb.edu/dokuwiki/doku.php?id=small_printer&rev=1639430782

Last update: 2021/12/13 21:26

